

Real-time 3D analytical FIB-SEM

**HITACHI**  
Inspire the Next

# NX9000

Real-time 3D analytical FIB-SEM



# FIB-SEM System for True 3D Structural Analysis

The analysis of internal structures are important for a comprehensive evaluation of advanced materials and devices. The combined Scanning Electron Microscope (SEM) and Focused Ion Beam (FIB) system enables site-specific cross sectioning and lamella preparation, which have become indispensable for SEM observation and Transmission Electron Microscope (TEM) analysis.

The newly developed FIB-SEM system from Hitachi, the NX9000 incorporates an optimized layout for true high resolution serial sectioning to tackle the latest demands in 3D structural analysis and for TEM analysis.

The NX9000 FIB-SEM system allows the highest precision in material processing for a wide range of areas relating to advanced materials, electronic devices, biological tissues and a multitude of other applications.



Hitachi patents: US6118122, US6538264, US6828466, US7139698, US7345289, US7397050, US7397051, US7442845, US7525188, US7660759, US8198803, US8500719, US8642958, US8664588, and other patents, as of June, 2015

# Real-time 3D analytical FIB-SEM **NX9000**

## Features

SEM column and FIB column are orthogonally arranged to optimize column layout for 3D structural analysis.

Combination of high brightness cold field emission electron source and high sensitivity optics support analysis of a wide range of materials from magnetic materials to biological tissues.

Micro-sampling system and Triple Beam system allow high quality sample preparation for TEM and atom probe applications.

## Applications

Ion milling and observation at normal incidence in real-time for true analytical imaging ..... P.3

Cut & See•3D-EDS <sup>(1)</sup>•3D-EBSD <sup>(2)</sup> available for a wide variety of materials ..... P.4

High resolution, high contrast SEM column  
by high brightness cold cathode field emission electron source and high sensitivity optics ..... P.5

High resolution 3D structural analysis enabled by highly stable,  
high precision sample stage ..... P.6

High quality TEM•atom probe sample preparation using Micro-sampling<sup>®</sup>  
and Triple Beam<sup>®</sup> system ..... P.6

(1) Energy Dispersive x-ray Spectroscopy, option

(2) Electron BackScatter Diffraction, option

# Ion milling and observation at normal incidence in real-time for true analytical imaging

The SEM column and FIB column are orthogonally arranged to realize normal incident SEM imaging of FIB cross sections. Orthogonal column arrangement eliminates aspect deformation, foreshortening of cross sectional images and shift of FOV (Field Of View) during serial section imaging, which cannot be avoided by conventional FIB-SEM systems. The NX9000 images produced enable highly accurate 3D structural analysis. Optical correlative microscopy can be applied easily due to the benefit of surface planar EM imaging.

**Conventional FIB-SEM**

Serial section images with deformation and FOV shift due to oblique section imaging

SEM incident angle: 54°  
Deformed aspect ratio due to oblique section imaging

SEM incident angle: 54°  
After aspect ratio correction

**Orthogonally arranged FIB-SEM**

Serial section images without deformation or FOV shift thanks to normal incident imaging

SEM incident angle: 90°  
No deformation thanks to normal incident SEM imaging

Sample: Mouse brain neuron  
Sample courtesy of Yoshiyuki Kubota, Ph.D.,  
Neural Information Processing Systems (NIPS)

# Real-time 3D analytical FIB-SEM **NX9000**


**Cut & See** · 3D-EDS\* · 3D-EBSD\* available for a wide variety of materials

## Cut & See

Cut & See supports high resolution, high contrast imaging of biological tissues, semiconductors, and magnetic materials such as steel and nickel at low accelerating voltage.

Serial section images can be collected with high throughput due to the proper geometry of the ion and electron column.

1 nm step Cut & See

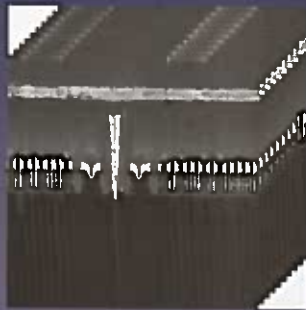

Serial section images



Reconstructed XZ plane



Sample : NAND flash memory  
SEM accelerating voltage : 1 kV  
Cutting interval : 1 nm  
Number of cut : 300




## 3D-EDS\*

Serial section SEM images and serial section elemental maps can be collected using 3D-EDS.

Large area silicon drift detector reduces acquisition time and enables elemental mapping at low accelerating voltage.


Sample : Fuel cell electrode  
SEM accelerating voltage : 5 kV  
Cutting interval : 100 nm  
Number of cut : 212

Sample courtesy of Prof. Naoki Shikazono, University of Tokyo

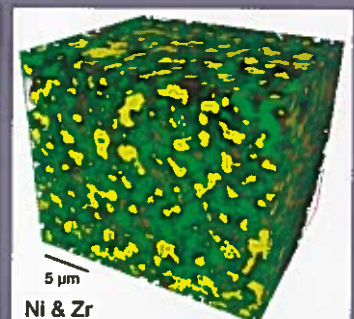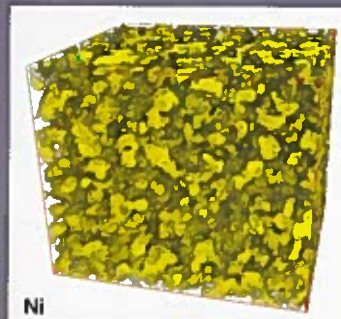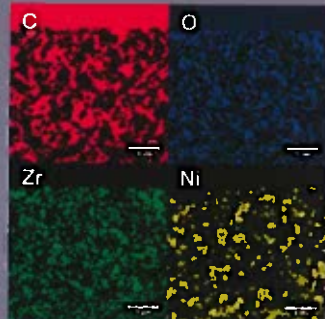


## 3D-EBSD\*

Simultaneous SEM, FIB and EBSD signals are obtained for 3D-EBSD without moving the stage during FIB sectioning and EBSD analysis. Accuracy and throughput of 3D crystal orientation analysis and segmentation yield high quality and less post-processing correction.

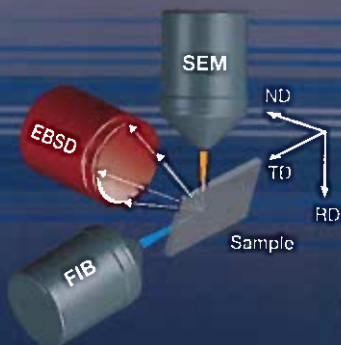


Sample : Ni  
SEM accelerating voltage : 20 kV  
Cutting interval : 150 nm  
Number of cut : 150

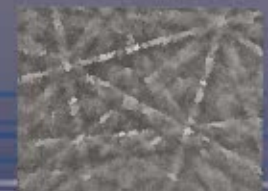

## High resolution, high contrast SEM column

by high brightness cold cathode field emission electron source and high sensitivity optics

Small source size and low energy spread of cold cathode field emission electron source allow high resolution imaging even at low accelerating voltage. In-column SED, In-column BSED, Retractable BSED\*, and STEM<sup>®</sup> detectors\* are placed at optimal positions to support a diverse range of samples and analytical purposes. Each signal can be simultaneously acquired at beam coincident point. For STEM imaging, bright field STEM detector\* and dark field STEM detector\* are available. FIB prepared sample for TEM and atom probe can be examined rapidly without taking sample out of the chamber.

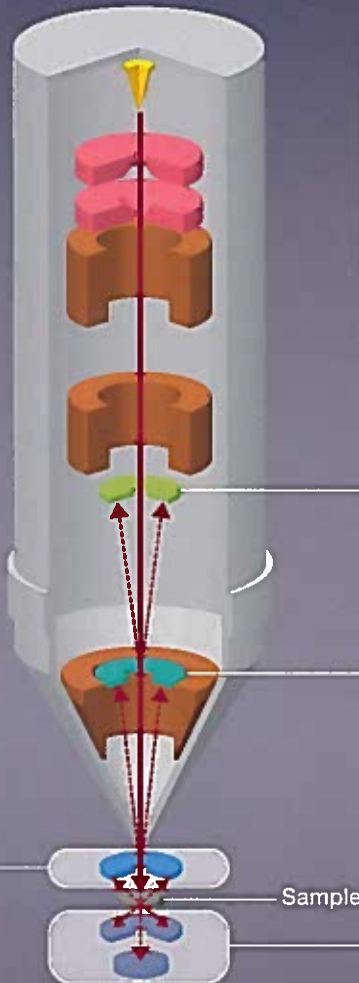

**In-column SED (Secondary Electron Detector)**

Fuel cell electrode

0.3 kV

0.1 kV

Ni ZrO Ni ZrO

**In-column BSED (Backscattered Electron Detector)**

Fuel cell electrode

0.3 kV

Mouse brain neuron tissue

1.5 kV

Ni ZrO

**Retractable BSED\***

Gallium nitride (GaN)

7 kV

**STEM detector\***

3D NAND flash memory

30 kV bright field

30kV dark field

# Real-time 3D analytical FIB-SEM NX9000


## High resolution 3D structural analysis enabled by highly stable, high precision sample stage

Optimized sample stage and sample holder significantly reduce drift for lengthy acquisition sessions. Together with stable ion source and optics, high precision serial FIB slicing is achieved to support high resolution 3D structural analysis at nanometer scale.


3D reconstruction



Sample : 3D NAND flash memory  
SEM accelerating voltage : 1 kV  
Cutting interval : 2 nm  
Number of cut : 498





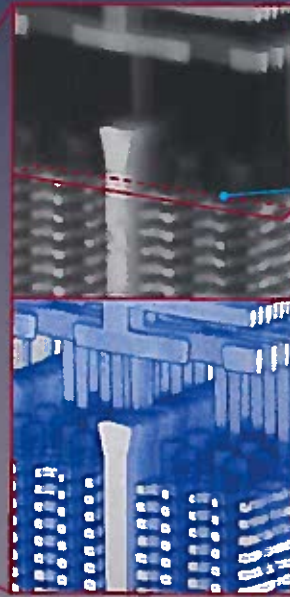


Poly Si layer measurement

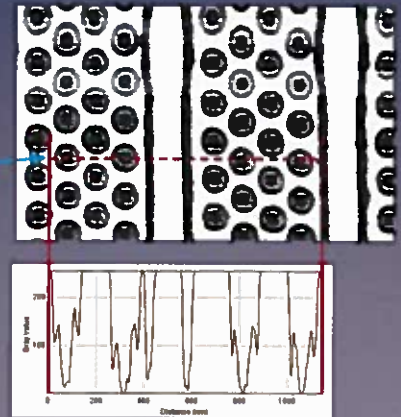

## High quality TEM·atom probe sample preparation using Micro-sampling\* and Triple Beam\* system

Micro-sampling system\* enables site-specific extraction of micro sample inside the chamber while Triple Beam system\* drastically reduces FIB-induced damage using Ar ion beam. For atom probe sample preparation, EBSD\* signal can be utilized for end point detection simultaneously. Needle shape sample can be prepared with specific grain boundary preserved which is difficult to identify by SEM.

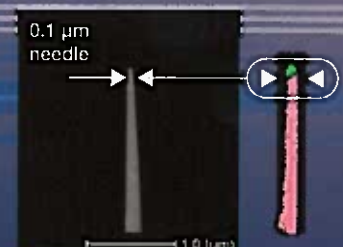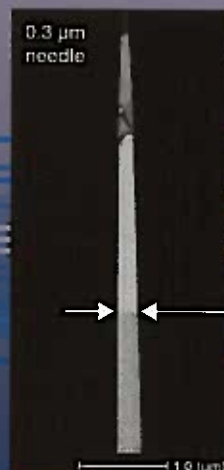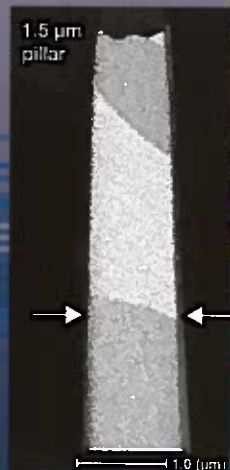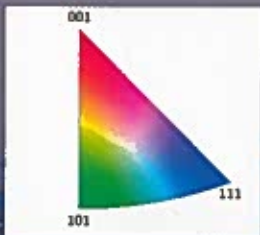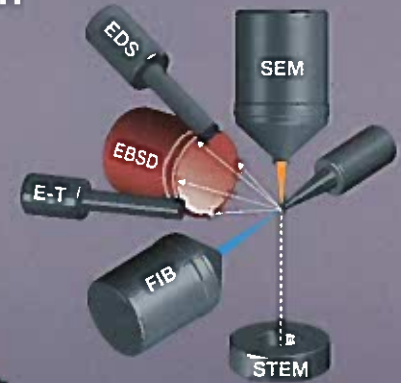

Sample : Metal material  
SEM accelerating voltage :  
1 kV for imaging  
20 kV for EBSD analysis

Option

## Specifications

NX9000		
SEM	Electron source	Cold cathode field emission source
	Accelerating voltage	0.1 - 30 kV
	Resolution	2.1 nm@1 kV 1.6 nm@15 kV
FIB	Ion source	Ga liquid metal ion source
	Accelerating voltage	0.5 - 30 kV
	Resolution	4.0 nm@30 kV
	Maximum probe current	100 nA
Standard detector		In-column SED / In-column BSED / Chamber SED
Stage	X	0 - 20 mm*1
	Y	0 - 20 mm*1
	Z	0 - 20 mm*1
	$\theta$	0 - 360°*1
	$\tau$	-25 - 45°*1
Maximum sample size		6 mm x 6 mm, 2 mm thick

\*1 Stroke is limited by each sample holder.

## Options

Ar ion 3 <sup>rd</sup> column	For low energy Ar ion beam polishing with real-time SEM end point detection functions.
Micro-sampling system	For micro sample extraction from bulk sample.
Multi gas injection system	For deposition of multiple gases. Gas needs to be specified.
•Carbon deposition gas	
•Tungsten deposition gas	
•Platinum deposition gas	
Retractable BSED	For low angle BSE detection. Retractable STEM detector cannot be mounted simultaneously.
STEM detector (STEM Type I)	For bright field STEM imaging.
Retractable STEM detector (STEM Type II)	For dark field STEM imaging. Retractable BSED cannot be mounted simultaneously.
Cold trap	Effective for reducing contamination.
Air protection sample transfer	Sample environment protection during transfer.
Side entry auto stage	TEM holder stage for direct transfer. Occupies EBSD port.
Air protection cooling holder	Side entry auto stage required.
Sample rotation holder	Side entry auto stage required.
Plasma cleaner	For reduction of hydrocarbon in SC chamber.
EDS	For X-ray Dispersive Spectroscopy.
EBSD	For crystal orientation analysis and mapping. Side entry auto stage cannot be mounted simultaneously.
A-TEM	For automatic TEM sample preparation.
Vector scan	For customised patterning by vector scan.
3D-EDS software	EDS system required.
3D-EBSD software	EBSD system required.
Chiller	For SEM lens system cooling.

NOTICE: For correct operation, follow the instruction manual when using the instrument.

Specifications in this catalog are subject to change with or without notice, as Hitachi High-Technologies Corporation continues to develop the latest technologies and products for our customers.

Copyright (C) Hitachi High-Technologies Corporation 2015 All rights reserved.

## Hitachi High-Technologies Corporation

Tokyo, Japan

[hitachi-hightech.com/global/science/](http://hitachi-hightech.com/global/science/)

24-14, Nishi-Shimbashi 1-chome, Minato-ku, Tokyo 105-8717, Japan

## Hitachi High-Tech Science Corporation


<http://www.hitachi-hightech.com/global/hhs/>

Head office

24-14, Nishi-Shimbashi 1-chome, Minato-ku, Tokyo 105-0003, Japan



## Layout

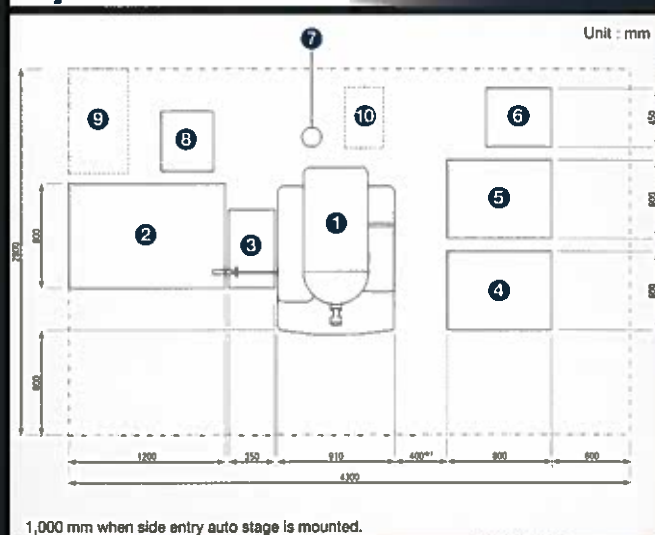

Unit	Width x length x height, mm	Weight, kg
① Main unit	910 x 1,280 x 2,015	1,500
② Operation table	1,200 x 800 x 720	60
③ Sub table	350 x 600 x 700	35
④ SEM control cabinet	600 x 800 x 2,000	350
⑤ FIB control cabinet	600 x 800 x 2,000	485
⑥ Transformer	450 x 500 x 450	70
⑦ Weight	Φ150 x 120	20
⑧ Scroll dry pump	500 x 460 x 319.5	35
⑨ Chiller (recommended model)*2	377 x 500 x 620	45
⑩ Argon cabinet**	290 x 465 x 823	50

\*2 Option

## Utilities

Room temperature	Set value	22°C±3°C
	Tolerance	Set value ±1°C
	Fluctuation	0.5°C/h or less
Humidity	Tolerance	35 - 60% no condensation
	Power	Single phase AC200/208/220/230 V±10% 30 A (50/60 Hz)
Grounding	D-class	100 Ω or less
	Cooling water	Flow
Pressure		50 - 100 kPa
Temperature		15 - 25°C
Temperature fluctuation		±0.5°C or less/10 min
N <sub>2</sub> gas	Purity	99.95% or higher
	Pressure	0.5 - 0.7 MPa